SOPORTE TÉCNICO

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For research use only. Not for human, veterinary or diagnostic use.

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TECHNICAL SUPPORT

Laboratory receipt, storage and reconstitution sheet

General procedure for CanLab México lyophilised reference materials. It complements — and does not replace — the laboratory's internal protocol, the lot COA and the compound SDS.

1. Receipt and inspection

- Match the lot number printed on the vial label against the certificate of analysis (COA) and the commercial invoice: all three must agree.
- Check the aluminium seal and stopper are intact; do not accept vials with a lifted seal or damaged glass.
- Record the reading of the packaging temperature monitor and keep it in the lot file.
- Notify any non-conformity in writing within 10 business days, quoting lot number and photographic evidence.

2. Storage

- Lyophilised: store at -20°C , protected from light and moisture, in the original closed vial.
- The powder is stable at ambient temperature in transit; long-term storage must be in a freezer.
- Before opening, let the closed vial equilibrate 15–20 minutes at room temperature to avoid condensation on the powder.

3. Reconstitution

- Solvent: bacteriostatic water or sterile reagent-grade water, according to the laboratory protocol. For compounds poorly soluble in water consult the COA and the compound sheet.
- Add the solvent slowly down the wall of the vial, without directing the stream onto the powder.
- Do not shake. Swirl the vial gently until the solution is clear.
- The working-solution concentration follows from the vial's net content and the solvent volume added (table).

WORKING-SOLUTION CONCENTRATION (MG/ML)

VIAL CONTENT	+ 1 ML	+ 2 ML	+ 3 ML
5 mg	5.00 mg/mL	2.50 mg/mL	1.67 mg/mL
10 mg	10.00 mg/mL	5.00 mg/mL	3.33 mg/mL

Concentration = vial net content ÷ solvent volume added. Nominal values; the actual content is the one declared on the lot COA.

4. Solution stability

- Keep the reconstituted solution at 2–8 °C, protected from light, and write the reconstitution date on the vial.
- Use within weeks, depending on compound and solvent; consult the compound sheet.
- Avoid freeze–thaw cycles. If the protocol requires it, aliquot into low-binding tubes and freeze once only.

5. Handling and protective equipment

- Nitrile gloves, safety glasses and lab coat; fume hood or cabinet when weighing powder.
- Do not ingest. Do not apply to people or animals. Laboratory use by qualified personnel only.
- Consult the compound SDS where available.

6. Waste

- Dispose of vials, solvents and contaminated material under local laboratory chemical-waste rules.
- Do not pour down the drain or mix with household waste.

7. Traceability

- Keep the COA, invoice and temperature-monitor record together with your laboratory's lot register.
- Note date of receipt, date of opening, solvent and volume used.

Solo para investigación. No apto para uso humano, veterinario ni diagnóstico.
For research use only. Not for human, veterinary or diagnostic use.

Every material in this catalogue is supplied exclusively for in-vitro and preclinical research in qualified laboratories. They are not drugs, are not approved by COFEPRIS, INVIMA, ANVISA or Health Canada for human or veterinary use, and must not be used in diagnostic procedures or administered to people or animals outside an authorised research protocol.